

# Ampere's Law Accessory

EM-6720



## Introduction

The Ampere's Law Accessory is used in combination with wireless rotary motion and magnetic field sensors and a current-carrying coil to verify Ampere's Law:

$$\oint \overrightarrow{B} \cdot \overrightarrow{dl} = \mu_o N I_{enclosed}$$

Detailed experiment instructions are available to download at <a href="https://www.pasco.com">www.pasco.com</a>. Type EX-5552 into the search to go to the Ampere's Law Experiment.

# **Equipment**

Ampere's Law Accessory EM-6720 includes:


1	Platform - Left
1	Platform - Right
1	Sensor Bracket
2	1/4-20 x 3/4" Thumbscrews
1	1/4-20 x 1/2" Brass Thumbscrew
6	6-32 x ¼" Brass Thumbscrews

Additional equipment required to perform the complete experiment (EX-5552):


1	Wireless 3-Axis Magnetic Field Sensor	PS-3221
1	Wireless Rotary Motion Sensor	PS-3220
1	Field Coil (500-Turn)	EM-6723
1	Zero Gauss Chamber	EM-8652
1	850 Universal Interface (or other DC power supply that outputs at least 1 A at 15 V	UI-5000
	PASCO Capstone Software	

# **Platform Assembly**


 Slide the coil into the slot in the right side of the platform.

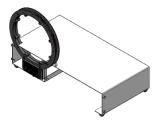


2. Secure the coil with the  $\frac{1}{4}$  x 20 brass screw that is stored in the platform.

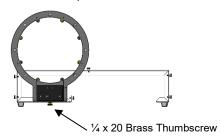

3. Remove the six 6-32 brass screws from the right side of the platform. Slide the left side of the platform into the right side, with the right side overlapping the left side.

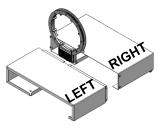


4. Secure the right side to the left side using the six 6-32 brass screws.

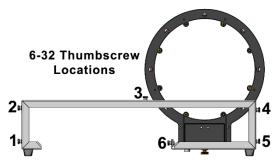

# **Sensor Assembly**

- Remove the two black thumbscrews from the Sensor Bracket.
- 2. Put the O-ring (included with PS-3220) on the largest step of the 3-step pulley and put the pulley on the Wireless Rotary Motion Sensor (PS-3220).
- Use one thumbscrew to attach the Wireless Rotary Motion Sensor to the Sensor Bracket.

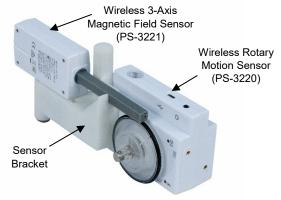

 Position the probe of the Wireless 3-Axis Magnetic Field Sensor (PS-3221) over the pulley of the Rotary Motion Sensor. Then secure it in place with the other black thumbscrew.

## **Technical Support**

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific 10101 Foothills Blvd. Roseville, CA 95747-7100

Phone: +1 916 462 8384 (worldwide)

8700-772-8700 (U.S.)

Web: www.pasco.com
Email: support@pasco.com

The Reference Guide will be updated periodically. For the latest revision of this Reference Guide, visit the PASCO Web site at

#### www.pasco.com/manuals

and enter the product number, EM-6720, in the text window.

## **Replacement Parts**

For information about possible replacement parts, contact Technical Support:

#### **Limited Warranty**

For a description of the product warranty, see the PASCO catalog. For more information visit www.pasco.com/legal.

#### Copyright

This PASCO scientific *Reference Guide* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited. (Revised 02/11/19)

#### Trademarks

PASCO, PASCO scientific, PASCO Capstone, and SPARKvue are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of their respective owners. For more information visit www.pasco.com/legal.

#### **Product End of Life Disposal Instructions:**

This electronic product is subject to disposal and recycling regulations that vary by country and region. It is your responsibility to recycle your electronic equipment per your local environmental laws and regulations to ensure that it will be recycled in a manner that protects human health and the environment. To find out where you can drop off your waste equipment for recycling, please contact your local waste recycle/disposal service, or the place where you purchased the product.

