
RADIATION CANS

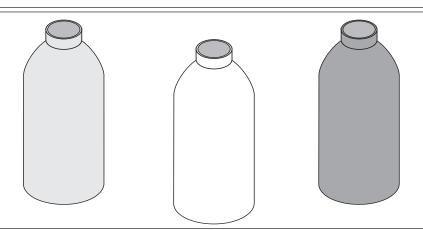

Introduction

The PASCO Radiation Cans are a set of three cans: one silver, one white, and one black. These cans are very useful in demonstrating that objects of the same material, but with different colors, will absorb and radiate heat at different rates. By filling the cans with water, and determining the temperature of each, one can confirm that the darker can heats and cools more quickly than the cans of lighter color.

Equipment

• Radiation Cans

Additional Equipment Recommended:

- Digital Thermometer, PASCO Model SE-9086 or, CI-6505A Temperature sensor with Science Workshop
- fan (optional)

Copyright Notice

The PASCO scientific 012-06658 instruction sheet is copyrighted and all rights reserved. However, permission is granted to non-profit educational institutions for reproduction of any part of the (Radiation Cans) instruction sheet providing the reproductions are used only for their laboratories and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Limited Warranty

PASCO scientific warrants the product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. The warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment after repair will be paid by PASCO scientific.

Address:PASCO scientific

10101 Foothills Blvd. P.O. Box 619011 Roseville, CA 95678-9011

Phone: (916)786-3800 FAX: (916)786-8905

email: techsupp@pasco.com web: www.pasco.com





Radiation Cans 012-06658A

Technical Support

Feedback

If you have any comments about the product or manual, please let us know. If you have any suggestions on alternate experiments or find a problem in the manual, please tell us. PASCO appreciates any customer feedback. Your input helps us evaluate and improve our product.

To Reach PASCO

For technical support, call us at 1-800-772-8700 (toll-free within the U.S.) or (916) 786-3800.

fax: (916)786-3292

e-mail: techsupp@pasco.com

web: www.pasco.com

Contacting Technical Support

Before you call the PASCO Technical Support staff, it would be helpful to prepare the following information:

If your problem is with the PASCO apparatus, note:

- Title and model number (usually listed on the label);
- Approximate age of apparatus;
- A detailed description of the problem/sequence of events (in case you can't call PASCO right away, you won't lose valuable data);
- If possible, have the apparatus within reach when calling to facilitate description of individual parts.

If your problem relates to the instruction manual, note:

- Part number and revision (listed by month and year on the front cover);
- Have the manual at hand to discuss your questions.

